

Getting start with Qsep₁



Packing List
包裝清單

C105200 (S1 & S2 Cartridge Kit)
C105200 (S1 & S2 卡匣組)

Qsep₁ power adaptor
Qsep₁ 電源供應線

Markers
標準品

C105200 (S1 & S2 Cartridge Kit)
C105200 (S1 & S2 卡匣組)



Qsep₁ power adaptor
Qsep₁ 電源供應線

Sample trays
(12 and 8 wells)
and Ethernet cable
樣品盤
(12和8孔)
及網路線

Qair_{box} and its
Power adaptor
Qair_{box} 及
電源供應線

Install Qsep₁ 安裝 Qsep₁

1. Unpack Qair_{box} and tighten the air tube with connector
將 Qair_{box} 開箱並把氣管插入轉接頭

Connector



2. Plug power cord into Qair_{box}
(* Please check the Qair_{box} label on the power cord)
請先將 Qair_{box} 接上電源
(* 請確認電源線上的標籤是屬於 Qair_{box})



label

3. Plug the power cord and Ethernet cable into Qsep₁
(Please make sure the label on the power cord is labeled Qsep₁)
將網路線及電源供應線接上 Qsep₁ (請確認電源線上的標籤是屬於 Qsep₁)



4. Connect the other side of Ethernet cable to the operation computer.
將網路線另一端連接您操作的電腦



5. Plug another side of air tube into Qsep₁ and turn on the Qair_{box}
連結氣管線並開啟 Qair_{box} 電源



Your Qsep₁ is READY TO USE 隨時可以使用您的 Qsep₁

***DO NOT switch on the instrument immediately after powering off
Wait at least 5 seconds**